

PRE TREATMENT OF SEDIMENT FOR GRAIN SIZE ANALYSIS WITH A LASER MICROGRANULOMETER (carbonates, organic matter, biogenic silica)

MATERIAL

1. Glassware:
Porcelain crucible, 50 ml Falcon® tube, glass slide, glass stick, wash bottle 100 ml. (volumetric flask for chemical preparations)
2. Equipment:
Laboratory weighing scale, centrifuge, bain-marie, vortex, microscope, laboratory oven, magnetic agitator.
3. Purified water: **H₂O_d**
4. Chemical:
 - Hydrochloric acid 37% [\[7647-01-0\]](#) : **HCl_A** 4% solution, **HCl_B** 20% solution
 - Hydrogen peroxyde 35% p/p (133vol.) [\[7722-84-1\]](#) : **H₂O₂** at 35% p/p (can be diluted if too brutal reaction)
 - Sodium carbonate [\[497-19-8\]](#) : **Na₂CO₃** 2M solution
 - Sodium hydroxyde [\[1310-73-2\]](#) : **NaOH** 1M solution
 - Sodium hexametaphosphate (calgon) [\[6891-31-1\]](#) : **calgon** 1g/L solution
 - * { ▪ Ethanol [\[64-17-5\]](#) : used if the carbonate removal or organic matter destruction are too brutal (one pressure of wash bottle will stop or mitigate the reaction)

Caution: All chemical reactions should be performed under the fume hood and following all security requirement (lab coat, glasses and gloves). Used specific bin for reaction supernatants.

METHOD

1. Sample preparation:
 - Dry the sample in the laboratory oven at maximum 40°C for 24 hours. Let it longer if the sample is not completely dried.
 - Take between 1 and 2 g of sediment. Residue of the different destructions need to be sufficient for the micro-granulometry analysis (40 mg for fine sediment, 150 mg for large sediment)
2. Carbonate removal:
 - Put the withdraw in the porcelain crucible. Flood with **HCl_A** (*). Should observed an effervescence corresponding to CO₂ emission.
 - Mix using the glass stick. Let the reaction continue for 15 min.
 - Add between 1 and 2 ml of **HCl_A**. Wait a few seconds. Repeat the action until the reaction is over (no emission).
 - To control the efficiency of the reaction, add between 1 and 2mL of **HCl_B**. If not, repeat the action.
 - When the reaction is completed, put the sample into a Flacon® tube. Add 50 mL of **H₂O_d**. Vortex to unsettle the sediment.

- Centrifuge 7 min at 3000 rpm. Remove the supernatant and resuspend the pellet in **H₂O_d**. Repeat this step 2 to 3 times.
- Dry in the oven at 40°C and weight when dry. *Comparison with weight before treatment will indicate the percentage of carbonate within the sample.*

3. Organic matter removal

- Realise the treatment directly in the tube with open lid (emission).
- Caution, the reaction can be really brutal. Add a few drops of **H₂O₂** to evaluate the reaction intensity beforehand (*).
- Add 20 mL of **H₂O₂**. Carefully mix to resuspend the sediment.
- Put the tube in the bain-marie at 65°C for 6 hours. Mix occasionally to resuspend the sediment. Add 1 to 2 ml of **H₂O₂** if needed.
- After 6 hours, take the tube out. Resuspend the sediment and add 1 to 2 mL of **H₂O₂**. Let the reaction occur at room temperature. Mix occasionally and add **H₂O₂** if needed. *It can take several days before the reaction is completed.*
- Once the reaction is completed, add 50ml of **H₂O_d**. Vortex to resuspend the sediment.
- Centrifuge 7 min at 3000 rpm. Remove the supernatant and resuspend the sediment in **H₂O_d**. Repeat this step 4 to 5 times.
- Dry in the oven at 40°C and weight when dry. *Comparison with weight before treatment will indicate the percentage of organic matter within the sample.*

4. Biogenic silica removal, 2 ways:

A. Sediment with low silica content

- a. Reaction within the seal tube using sodium carbonate.
- b. Add 40 ml of **Na₂CO₃** (2M). Mix and put the tube in the bain-marie at 90°C for 6 hours. Mix every 2 hours.
- c. After 6 hours, add 50 ml of **H₂O_d**. Vortex and resuspend the sediment.
- d. Centrifuge 7 min at 3000 rpm. Remove the supernatant and resuspend in **H₂O_d**. Repeat this step 5 to 6 times. *Caution, ample formation of salt. Rinse abundantly.*
- e. Check on a smear slide at x500.
 1. If absence of diatom, dry in the oven at 40°C and weight when dry. *Comparison with weight before treatment will indicate the percentage of biogenic silica within the sample.*
 2. If diatoms, repeat previous step **b** to **e** with only 2 hours bain-marie. Dry in the oven at 40°C and weight when dry. *Comparison with weight before treatment will indicate the percentage of biogenic silica within the sample.*

B. Sediment with high silica content

- a. Reaction within a seal tube using sodium hydroxide 1M.
- b. Add 20 ml of **NaOH** (1M). Mix and put in the bain-marie at 70°C for 8 hours. Mix every 2 hours.
- c. After 8hours, add some **NaOH** and put back in the bain-marie at 70°C for again 8 hours. This step can be done several times.
- d. Once the reaction is completed, add 50 ml of **H₂O_d**. Vortex and resuspend the sediment.
- e. Centrifuge 7 min at 3000 rpm. Remove the supernatant and resuspend in **H₂O_d**. Repeat this step 5 to 6 times.
- f. Check on a smear slide at x500. No diatoms should be observed.
- g. Dry in the oven at 40°C and weight when dry. *Comparison with weight before treatment will indicate the percentage of biogenic silica within the sample.*

5. If the sediment is agglomerate, add 1 ml of **calgon** (1 g/L solution). Mix and let the reaction for 1 to 24 hours depending on the sample nature. It should be fully dissociated.
6. The sample is ready for grain size analysis with a laser microgranulometer.
MALVERN: MASTERSIZER hydro2000G with automatic sampler (0,020 to 2000µm).
Measurement according to Mie's theory with Fraunhofer's approximation. Results in percentages.

Literature :

Aufrère L. (1937). Introduction à l'Étude industrielle de la Silice en archéologie primitive, bulletin de la Société préhistorique de France, tome 34, n°3, 160-168

Borisovna Albot O. (2016). Pleistocene cyclostratigraphy on the continental rise and abyssal plain of the western Ross Sea, Antarctica, Master thesis Victoria University of, New Zealand

Lewis W.D., McConchie D. (1994). Analytical Sedimentology, Springer-Science + Business Media, B.V.(ed) 197,chapter 5, 61-84

Dron R. (1998). Mécanisme de la réaction alcali-silice, Bulletin des laboratoires des Ponts et Chaussées, 214, 61-68

Francus P., von Suchodoletz H., Dietze M.,V.Donner R., Bouchard F., Roy A.J., Fagot M., Verschuren D., Kröpelin S. (2013).Varved sediments of Lake Yoa (Ounianga Kebir, Chad) reveal progressive drying of the Sahara during the last 6100 years. Sedimentology Volume 60, 911-934

Passchier S. (2011). Linkages between East Antarctic Ice Sheet extent and Southern Ocean temperatures based on a Pliocene high-resolution record of ice-rafted debris off Prydz Bay, East Antarctica, Paleoceanography, vol. 26, PA4204